## Vanderbilt University Medical Center Research Laboratory Eyewash Log

Department	Location
-	

## Instructions:

- 1. All eyewash stations should be identified with a highly visible sign and should be free of debris.
- 2. Activate eyewash at least monthly to flush line and verify operation. Allow water to run for at least 30 seconds. Water should flow from both nozzles at equal velocity. There should be sufficient water pressure to deliver water at least 1.5" distance from the eyewash hardware.
- 3. If eyewash is not fully operational, call Facilities Management at 615-322-2041 or other applicable maintenance staff.

Date	Initials of staff performing test.	Easy to activate. (Y/N)	Delivers clear, tepid water. (Y/N)	Both sprays deliver equal velocity. (Y/N)	Stays on hands- free until turned off. (Y/N)	Nozzles are clean & covered when eyewash is not in use. (Y/N)